

Result 3



PLANETARY MONO MILL PULVERISETTE 6 CLASSIC LINE

main disk speed: 650 rpm

250 ml grinding bowl made of stainless steel

+ 15x20 mm grinding balls made of stainless steel.

Feed quantity: 70 g

Feed Size: < 5 mm

Grinding time: 8 min

Final fineness: fine

Comments: About 70g of cold cocoa beans fit into the 250ml bowl.

Within the first 3 minutes of grinding, the sample starts melting and the volume of sample collapsed.

After 8 minutes of grinding, no coarse particles are detectable between the fingers (so all particles use to be < 100µm).

The temperature of the sample is about 70°C.

A longer grinding is still possible.